



1321 S VALLEY CENTER DRIVE
BAY CITY, MI 48706
(989) 686-0350 FAX# (989) 686-0399
1-800-248-5060

Troubleshooting Steps for Goggle Sterilization Cabinets

1. How old is the cabinet (when was it purchased)?
2. What is the lot# off the bottom of the cabinet?
3. Does the bulb need replacing?
4. Check the power supply.
5. Make sure the timer is turned past 2.
6. Are both doors securely shut and locked
7. Is the bulb properly installed? Make sure both prongs on the bulb are in the light socket correctly and then twist until you feel a click
8. Make sure the starter is installed correctly. Twist and Click.

NOTE: DO NOT USE A 3 WIRE ADAPTER ON THE POWER CORD ON THIS UNIT!

WARNING: NEVER LOOK AT AN OPERATING ULTRAVIOLET LAMP WITH UNPROTECTED EYES!